



PRODUCT INFORMATION

MuA Transposase, conc.

Pub. No. MAN0013390

Rev. Date 24 August 2018 (Rev. B.00)

#F-750C

Lot _ Expiry Date _

Store at -25 °C to -15 °C

Components	#F-750C
MuA Transposase, conc., 1.1 µg/µL	22 µg

For Research Use Only. Not for use in diagnostic procedures.

www.thermofisher.com

Description

MuA Transposase is a single purified polypeptide that catalyzes transposition reaction *in vitro*.

Source

An *E.coli* strain that carries the cloned MuA gene from bacteriophage Mu.

Storage buffer

25 mM HEPES, pH 7.8, 0.1 mM EDTA, 2 mM DTT, 500 mM KCl, 50 % glycerol, non-ionic detergent.

Assay buffer for exonuclease and endonuclease reactions

25 mM Tris-HCl, pH 8.0 at 25°C, 5 mM MgCl₂, 5 mM NaCl.

Concentration was determined spectrophotometrically using Pierce 660 nm Protein Assay and Bovine Serum Albumin standard curve (0.125-0.5 mg/mL).

Notes

- Enzyme is supplied with a tube of Dilution Buffer (#F-776).
- It is not recommended to store enzyme dilutions prepared in the buffer.

Rev.5

Activity was determined by MuA Transposition Reaction/ Transformation Assay

Transposition reactions (20 µL) were performed using 1 ng/µL of the Entranceposon (M1-CamR), 18.5 ng/µL of the Control Target DNA (9243 bp, AmpR) and 22 ng/µL of the MuA Transposase in 1X reaction buffer (25 mM Tris-HCl pH 8.0 at 20 °C; 10 mM MgCl₂; 110 mM NaCl; 0.05 % Triton® X-100; 10 % glycerol). The reaction mixtures were incubated for 1 h at 30 °C followed by heat-inactivation of the MuA Transposase for 10 min at 75 °C. Transposition reaction volume of 10 µL was transformed into chemically competent *E.coli* cells using standard protocol (transformation efficiency > 10⁶ cfu/µg pUC57). Dilutions of the transformation mixture were plated on LB plates supplemented with 50 µg/mL ampicillin and 10 µg/mL chloramphenicol. As a result more than thousand chloramphenicol resistant colonies were recovered per single transposition reaction.

Exonuclease activity determined by Labeled Oligonucleotide (LO) Assay: No detectable degradation after incubation of single-stranded or double-stranded radiolabeled oligonucleotides with MuA Transposase.

Endonuclease Contamination: No detectable degradation was observed after incubation of supercoiled plasmid DNA with MuA Transposase.

Technical support

US: ts.molbio@thermofisher.com

Europe, Asia, Rest of World: ts.molbio.eu@thermofisher.com

Web: www.thermoscientific.com

Important Licensing Information:

This product may be covered by one or more Limited Use Label Licenses. By use of this product, you accept the terms and conditions of all applicable Limited Use Label Licenses.

PRODUCT USE LIMITATION

This product is developed, designed and sold exclusively *for research purposes and in vitro use only*. The product was not tested for use in diagnostics or for drug development, nor is it suitable for administration to humans or animals.

Please refer to www.thermofisher.com for Material Safety Data Sheet of the product.

© 2018 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries.